


Try Out

When the bot successfully boots, the images from the RGBD camera will show up on the screen. And all leds on the arm will turn green.

PLEASE make sure you WAIT for the leds to turn green. Do NOT move the arms before it's ready or people might get hurt.

Operate the arm

The robot is equipped with a mechanical arm base, allowing users to conveniently place the mechanical arm on the base to prevent it from moving randomly.

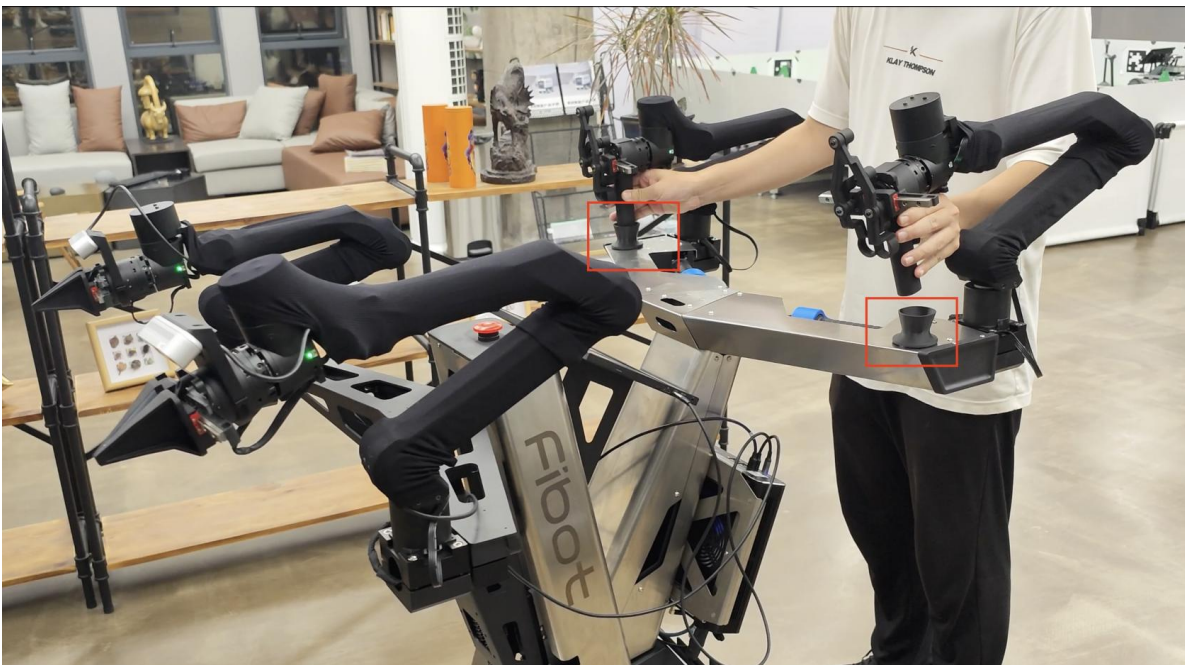

Move the robot around

To move the robot around, you don't have to use a remote(You can if you want though). Just grab the hand rail, and push/pull the robot anywhere you want. In free movement mode, you can move the robot forward and backward, left and right, and rotate it freely.



Use the arm lifter

We have designed a lifting platform for the mechanical arm to expand its workspace. You can use the blue rotary button on the handle to control the height of the lifting platform.

The lifting platform dramatically increased the workspace of the mechanical arm, enabling vertical opreation.





00 #5

0 Yangfan He 00 3 00 2024 08:38:23

0 Yangfan He 00 23 00 2024 13:44:02